

How to Use Disinfectant Cloth Wipes

This protocol is for the use of the reusable disinfectant wipes made by Colorado State University.

1. Remove a single cloth from the jar and make sure this cloth is wet/damp from the ethanol in the jar.
2. Wipe the desired surface with the wipe.
3. Allow the surface, that was just wiped, to air dry.
4. Put the cloth back into the jar unless it is grimy, at which point discard that wipe.

How to Use Disinfectant Spray Bottles

This protocol is for the use of 1% hydrogen peroxide (H₂O₂) spray bottles. This spray only requires minimal contact time, so the surface can be wiped with a paper towel after it is sprayed.

1. Spray the desired surface until it appears slightly damp.
2. Using a disposable paper towel, wipe down the surface.
3. Dispose of paper towel.